IDEXX **Quanti-Tray**

Insert and Most Probable Number (MPN) Table

Quanti-Tray Certificate of Sterility

This certifies that the enclosed Quanti-Trays have been sterilized with ethylene oxide.

For further information or documentation, contact IDEXX Laboratories, Inc.

North/South America: 1-207-556-4496/1-800-321-0207

Europe: 00800 4339 9111

UK: +44 (0) 1638 676800

China: +86-21-61279528

Japan: +81 422 71 5921

Australia: 1800 655 978

IDEXX

IDEXX Laboratories, Inc.
One IDEXX Drive
Westbrook, Maine 04092 USA

IDEXX Quanti-Tray

English Version

Introduction




IDEXX Quanti-Trays are designed to give quantitated bacterial counts of 100 mL samples using IDEXX Defined Substrate Technology reagent products. Add the reagent/sample mixture to a Quanti-Tray, seal it in a Quanti-Tray Sealer and incubate per the reagent instructions. Then count the number of positive wells and use the Most Probable Number (MPN) Table attached to determine the MPN.

Contents


This package contains sterile 51-well Quanti-Trays.

User Instructions

1. Use one hand to hold a Quanti-Tray upright with the well side facing the palm.
2. Squeeze the upper part of the Quanti-Tray so that the Quanti-Tray bends toward the palm.
3. Open the Quanti-Tray by pulling the foil tab away from the well side. Avoid touching the inside of the foil or tray.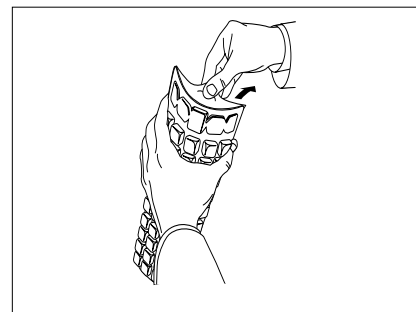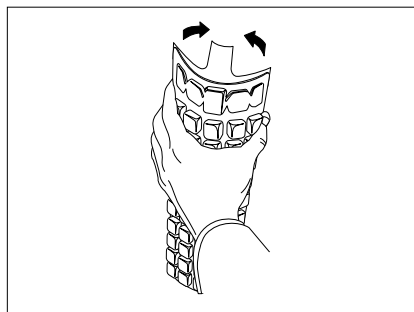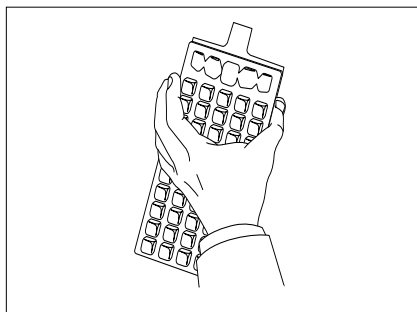

4. Pour the reagent/sample mixture directly into the Quanti-Tray, avoiding contact with the foil tab. Allow foam to settle.
5. Place the sample-filled Quanti-Tray onto the rubber tray carrier of the Quanti-Tray Sealer with the well side (plastic) of the Quanti-Tray facing down to fit into the carrier.
6. Seal according to the Quanti-Tray Sealer instructions.
7. Incubate according to reagent instructions.
8. Count positive wells and refer to the MPN table on the back of this instruction sheet to find the MPN.
9. Dispose of media in accordance with good laboratory practices.

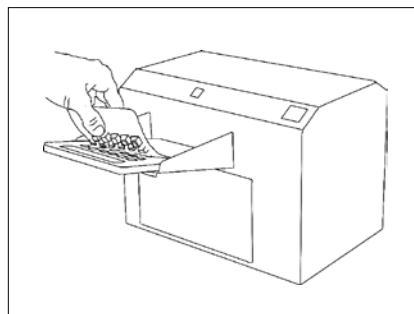For Technical Assistance
U.S./Canada 1-800-321-0207 or 1-207-556-4496
www.idexx.com/water

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